

Ulvac-PHI Versaprobe II XPS

The core technology of the VersaProbe II is PHI's patented, monochromatic, micro-focused, scanning x-ray source which provides excellent large area and superior micro-area spectroscopy performance. Standard features include PHI's patented dual beam charge neutralization technology, a floating column Ar ion gun for XPS sputter depth profiling, angle resolved XPS and a cold stage for temperature dependent studies.

Prerequisites for operating the Versaprobe tool:

- a) Obtain a NRF ID (if you do not already have one)
- b) Receive "one on one" training and certification from NRF Staff. Discuss your process with a staff member.

Safety



- **High Voltage** - High Voltage is used throughout the system. System maintenance may only be performed by NRF Staff. Do not remove any tool covers or defeat any interlock on this system.

1.0 Pre-Operation

1.1 None

2.0 Software Navigation

2.1 Start the Watcher and Smartsoft software if not already open

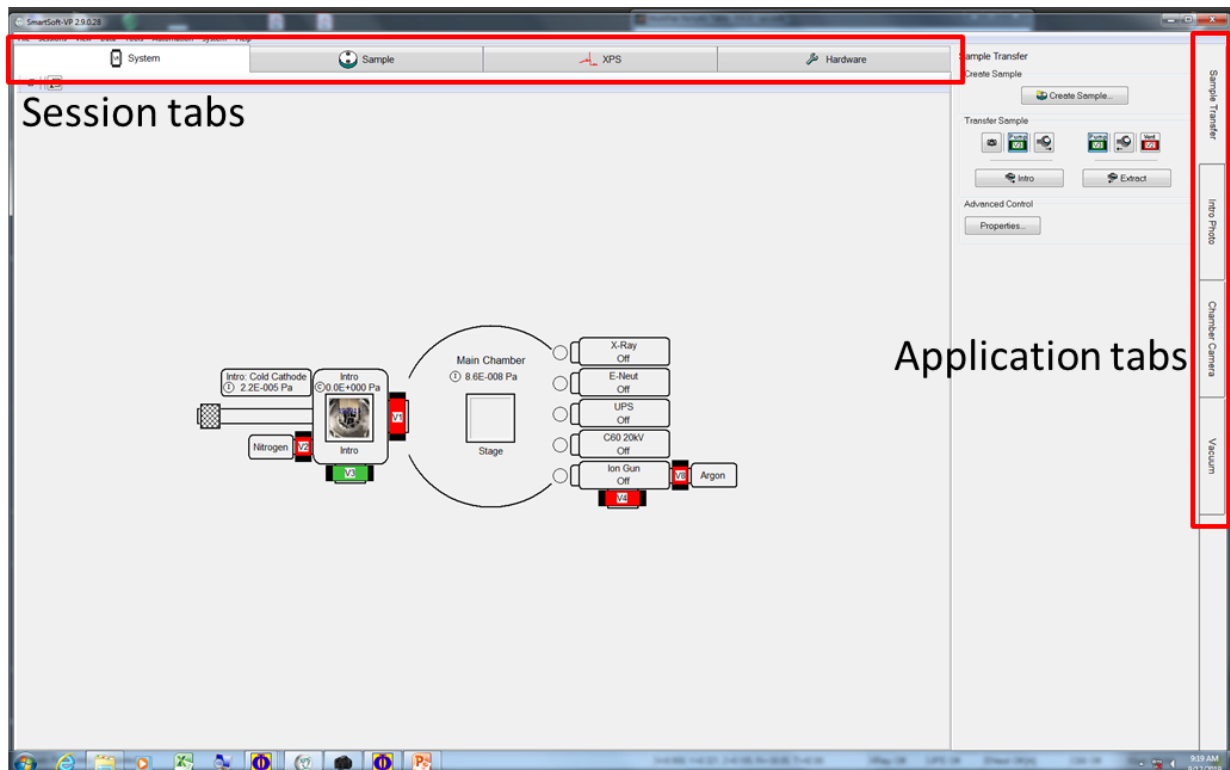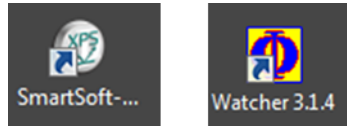


3.0 Sample Mounting

3.1 Samples must be mounted on a holder using the double-sided tape that is provided, or use the metal masks. Users can only use the 1 or 2 inch holders.

3.1.1. 1 inch holder - Normal use


3.1.2. 2 inch holder - For large samples or to analyze a lot of samples at once (*Stage tilt angle is fixed to 45 degrees when in use*)

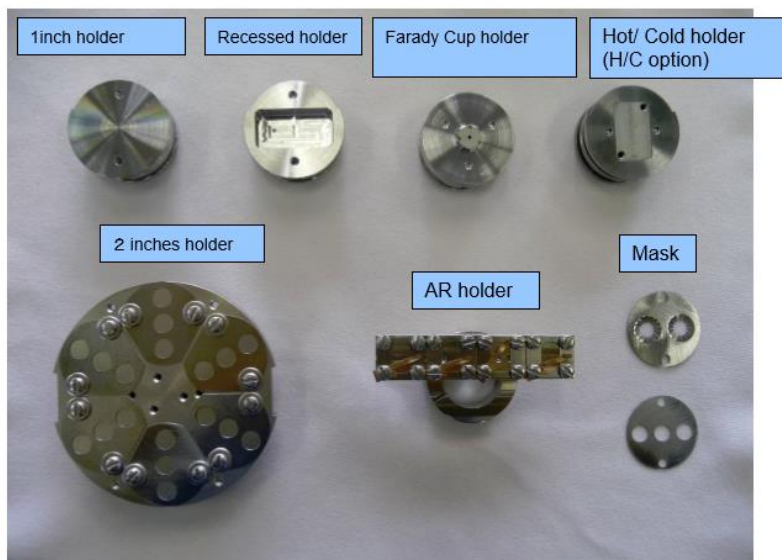


3.1.3. Do not touch your sample, the holder, or the business end of any tweezers with your bare fingers. Always use gloves and tweezers to mount and manipulate the sample and holder.

4.0 Sample Loading

- 4.1 Go to the 'System' session tab and click on the 'Sample transfer' application tab.
- 4.2 Click the 'Create sample' button and input a new platen name. The typical convention used is a YYMMDD format.
- 4.3 The 'Data Manager Properties' window will pop up and you can select the correct holder type (25mm, 60mm or Angle) from the pull-down menu.

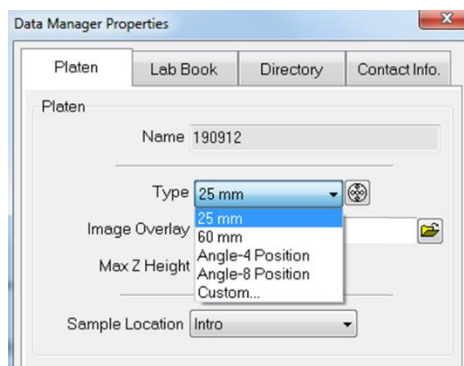


- 4.4 Click on the 'Directory' tab in the properties window to select, or create the location to store your data. All data must be stored in a subdirectory of C:\Data Files\Users. Use the 'Browse' button to open up a file explorer window and close the 'Properties' window when done.

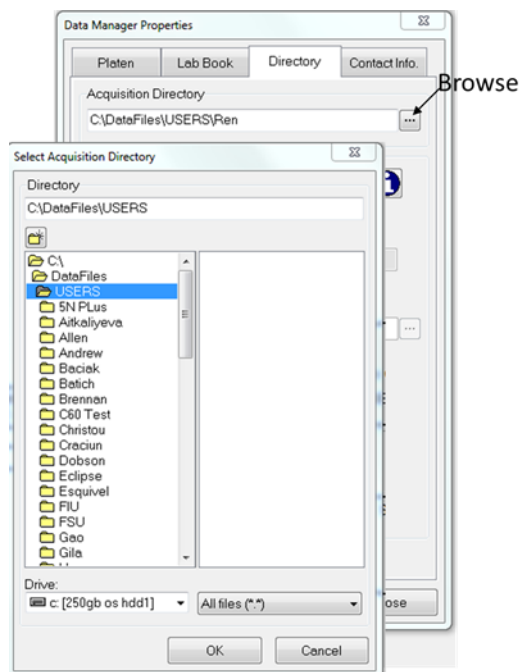


4.5 While still in the 'System' session tab, click on the 'Vacuum' application tab.

4.6 Select 'Backfill intro' from the 'Task' pull-down menu. Then click on 'Process Task'. This will vent the intro chamber (loadlock chamber).

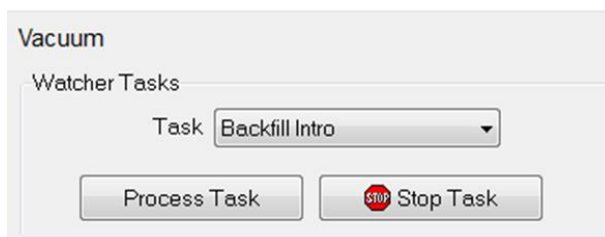


4.7 Remove the cap and place upside down on a surface. Do not put the sealing surface of the cap in contact with another surface.

4.8 Use the large tweezers hanging from the transfer arm to grab the holder and place it on the fork of the intro rod as shown in step 3 in the image below.

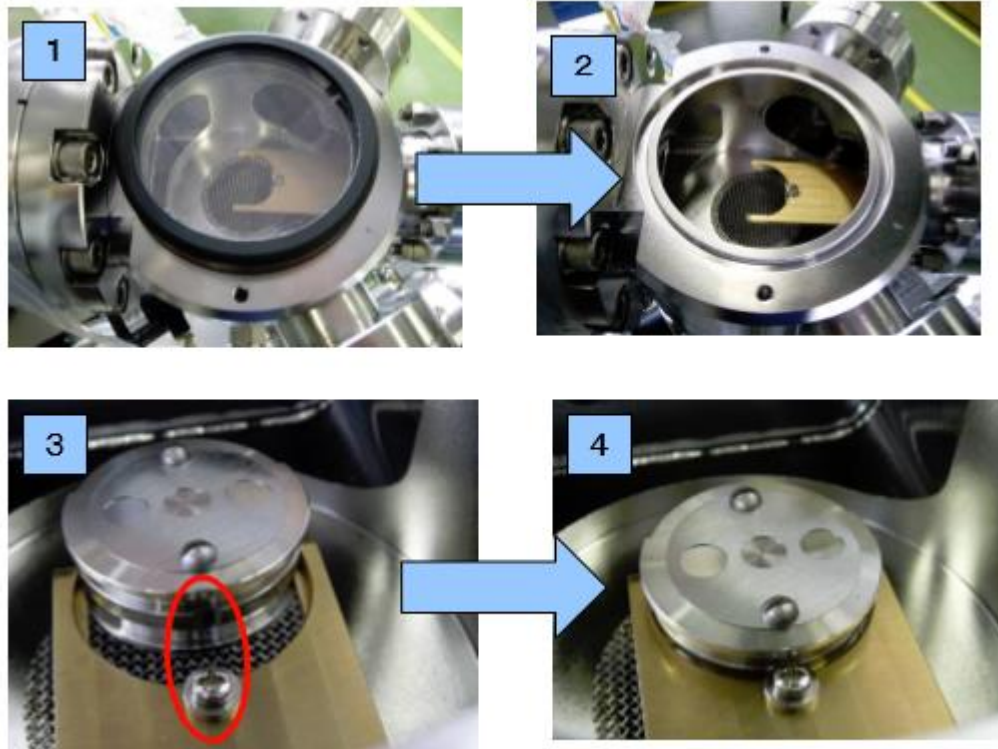


4.9 Ensure that the indented part of the holder is aligned to the screw as shown in the red circle. Slide the holder all the way back to the aligning screw as shown in step 4 above. Ensure that it is locked by trying to rotate the holder.


4.10 For the 2 inch holder, look through the side viewport and confirm that the arm forks are in the slot of the holder (like image A) and not just sitting on top (like image B). Also ensure that it is locked on the aligning screw by trying to rotate the holder.

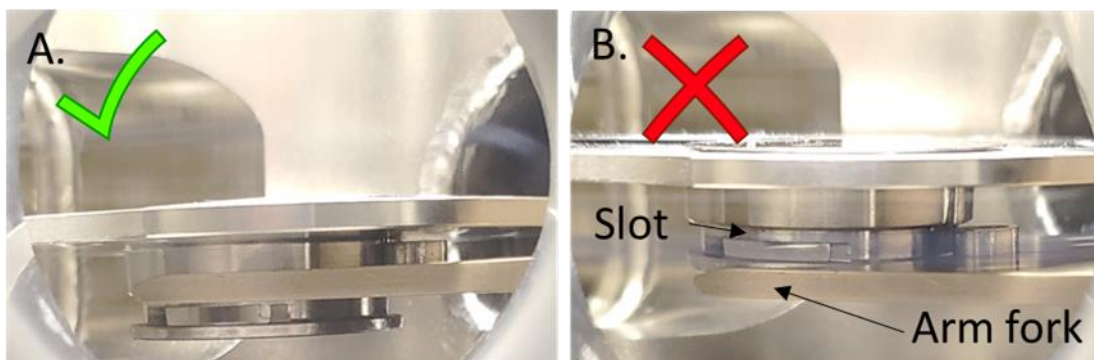

4.11 Leave the cover off and click the 'Intro' button.

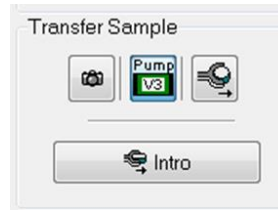

- 4.12 Click OK for the dialogue box. The Intro camera will take a photo.
- 4.13 Put the Intro lid back on, then click OK in the dialog box to start pumping.
- 4.14 After the specified pumping time, the stage moves to the transfer position and V1 opens then the dialog box below will open.
- 4.14.1 Watch inside the analysis chamber through the view port while slowly pushing the Transfer rod (red box) into the chamber until **you hit the stop collar** (red circle).
- 4.14.2 You can now click OK and the stage will move up to dock with the sample.

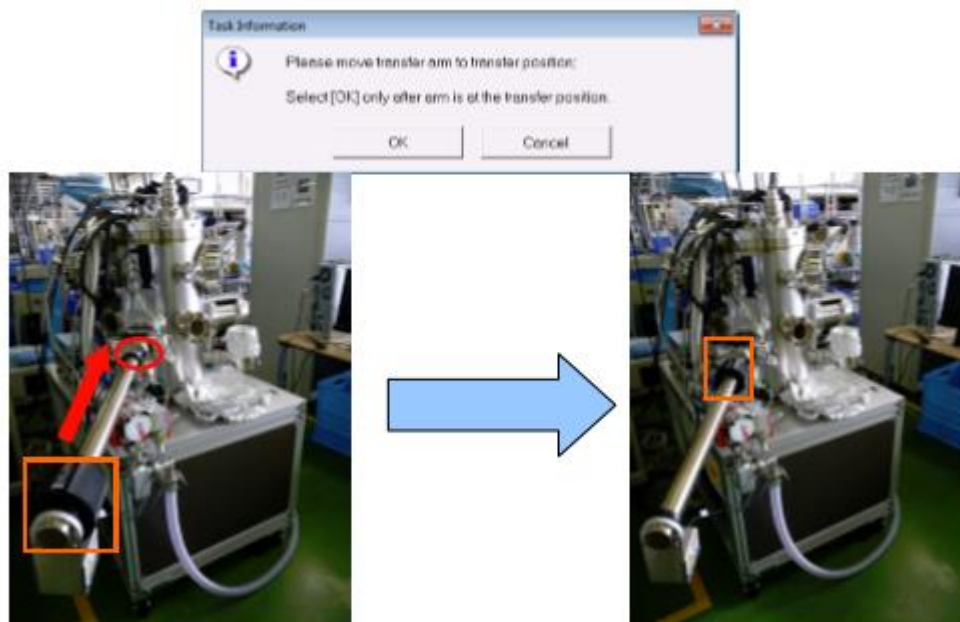


- 4.15 Watching through the viewing port, slowly pull the intro rod back completely. V1 should close automatically when the rod is completely withdrawn. Click the OK button in the message dialog to center the stage.
- 4.15.1 If you see that the sample is moving back with the arm, then push it all the way back in and try pulling out again.

5.0 Sample Set-up

5.1 Go to the 'Sample' session tab and click on the 'Z-align' application tab.

5.2 In the left region of the screen, you will see the photograph of the sample holder taken during the intro step. Place the cursor on a location you would like to analyze and right click on the mouse. Select 'Create point' from the menu.

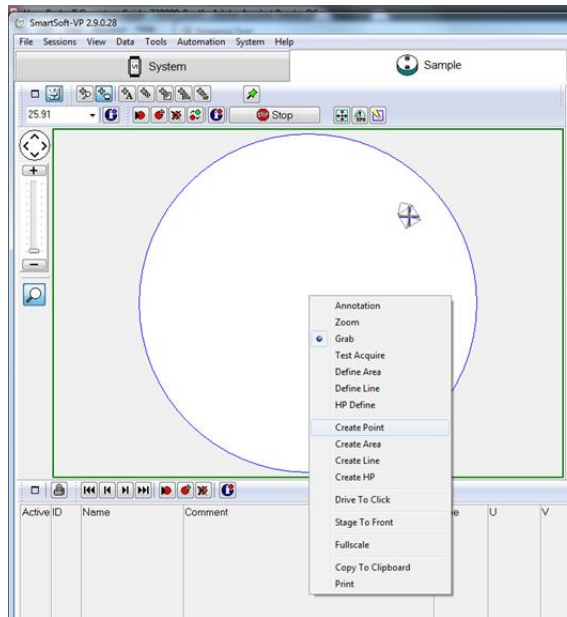

5.3 Measurement locations will be added to the table at the bottom. Repeat the process for other locations.

5.4 Z-alignment

5.4.1. From the position list, change the Z-position to 15mm for all locations and press the 'Update position' button for each one.

Drive to position Update position


Active	ID	Name	Comment	Type	U	V	Z	Rotation	Tilt
<input checked="" type="checkbox"/>	1	Sample 1		Point	-0.905	0.931	15.000	0.00	45.00
<input type="checkbox"/>	2	Sample 2		Point	5.374	-1.609	15.000	0.00	45.00

5.4.2. De-select all of the check marks in the 'Active' column on the list except for one.


5.4.3. Press the 'Drive to position' button

5.4.4. Press the 'Properties' button near the Z-align tab and graph window will appear.

5.4.5. Press the 'Z-align' button.

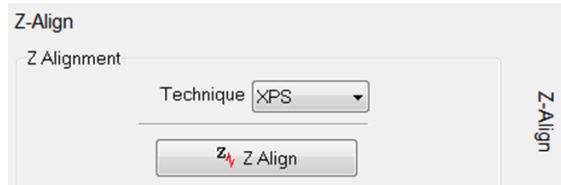

5.4.6. Ensure that the alignment was able to find the peak (green line passes through the peak) as shown below. It sometimes fails, so just press the 'Z-align' button again if it does.

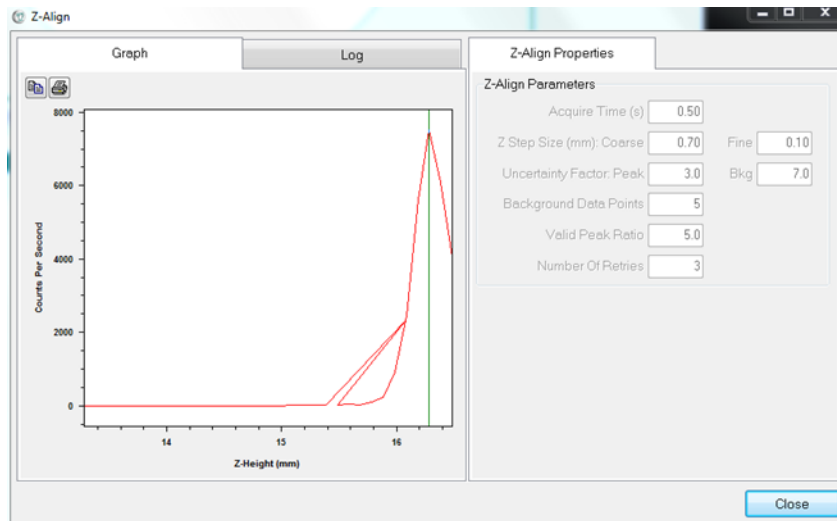

5.4.7. Press the 'Update position' button to save the new aligned position.

5.4.8. Repeat for all other locations.

6.0 XPS Spectrum

6.1 Go to the 'XPS' session tab and then the 'Spectrum' application tab.

6.2 Choose the Xray Setting from the drop-down menu that fits your analysis area and application

6.3 There are two modes to be used for analysis, Survey or Multiplex.

6.3.1. Survey - used for surface qualitative analysis (i.e. what elements are present)

6.3.2. Multiplex – used for quantitative or chemical state analysis

6.4 There are two ways to set up the survey and multiplex analysis regions.

- I. Go to the “regions” area of the screen and click on the periodic table
 - a. On the new periodic table, click the ‘Acquisition setup’ tab.
 - b. Select “Su” for a survey and select any elements that you wish to analyze
- II. Go to the “regions” area of the screen and use the plus/minus buttons to add or subtract region cells
 - a. Select the first cell and right click on the mouse, hover over “All elements”, and select “Su” for a survey
 - b. Do the same for all cells but select an element instead of Su

Spectrum

Spectrum Settings

TEMP

Load Save Delete File...

Spectrum

Spe More Add Q

Acquisition Directory

C:\DataFiles\USERS\Hays\delete

Source

Source FXS

Source Setting

X-Ray Setting 100u25W15KV

Spectrum Parameters

Analyzer Mode Scanned

Pass Energy 23.500

Time Per Step (ms) 20

Cycles 1

Analysis Time (min) 10.1 Runtime 11.2

Regions

+ -

Active Name Sweep Pass Lower Range

Active	Name	Sweep	Pass	Lower	Range
<input checked="" type="checkbox"/>	Su1s	10	93.900	0.0	1200.0
<input type="checkbox"/>	C1s	10	23.500	278.0	20.0
<input type="checkbox"/>	O1s	5	23.500	525.0	15.0
<input type="checkbox"/>	Si2p	10	23.500	94.0	20.0
<input type="checkbox"/>	N1s	10	23.500	391.0	20.0

Advanced Controls

XPS Setup...

Periodic Table

Acquisition Setup

Peak ID Filter

Acquisition Setup

Su 2 3 4 5 6 7 8

H Li Be B C N O F Ne

Na Mg Al Si P S Cl Ar

K Ca Sc Ti V Cr Mn Fe Co Ni Cu Zn Ga Ge As Se Br Kr

Rb Sr Y Zr Nb Mo Tc Ru Rh Pd Ag Cd In Sn Sb Te I Xe

Cs Ba La Hf Ta W Re Os Ir Pt Au Hg Tl Pb Bi Po At Rn

Fr Ra Ac

Ce Pr Nd Pm Sm Eu Gd Tb Dy Ho Er Tm Yb Lu

Th Pa U Np Pu Am Cm Bk Cf Es Fm Md No Lr

TEMP

Load Save Delete

Spectrum

Spe M

Acquisition Directory

C:\DataFiles\USERS\Hays\delete

Source

Source FXS

Source Setting

X-Ray Setting 100u25W15KV

Spectrum Parameters

Analyzer Mode Scanned

Pass Energy 23.500

Time Per Step (ms) 20

Cycles 1

Mn2p-641

Mn2p1-652

Mn2p3-641

Mn3p-48

Mn3s-83

Mn LMM-900

Mn LMM1-852

Mn LMM2-944

Regions

+ -


Active Name Sweep Pass Lower Range

Active	Name	Sweep	Pass	Lower	Range
<input checked="" type="checkbox"/>	Su1s	10	93.900	0.0	1200.0
<input type="checkbox"/>	C1s	10	23.500	278.0	20.0
<input type="checkbox"/>	O1s	5	23.500	525.0	15.0
<input type="checkbox"/>	Si2p	10	23.500	94.0	20.0
<input type="checkbox"/>	N1s	10	23.500	391.0	20.0

Advanced Controls

XPS Setup...

6.5 You can change the transition to be measured by right clicking on the element and choosing another transition. For example, you can choose to measure the Si2s or Si2p peaks in the figure below.

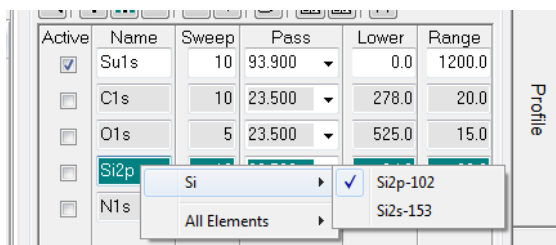

6.6 Set the analysis conditions.

6.6.1. Sweeps – The number of times to scan the energy range of interest. Typically 5-10

6.6.2. Set the energy range using the “lower, range, and upper” cells

6.6.3. Pass energy – Typically 93 for survey and 23.5 for multiplex. A higher number provides good signal to noise for qualitative analysis and lower numbers provide more accurate peak position for chemical state analysis.


6.6.4. Typically keep the defaults for ev/step and time per step

6.7 Click on the box under “Active” for the survey


6.8 Ensure that charge neutralization is set to automatic.

6.8.1. Click on the ‘XPS setup’ button.

6.8.2. Click on “Auto” for both E-neutralization and Ion gun.

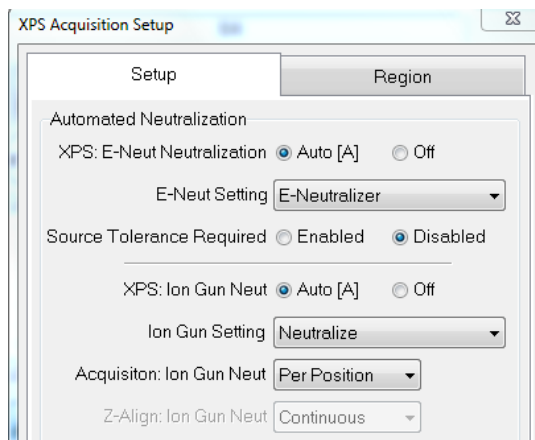

6.9 A pre-sputter can also be set up in this window to clean of any surface contamination.

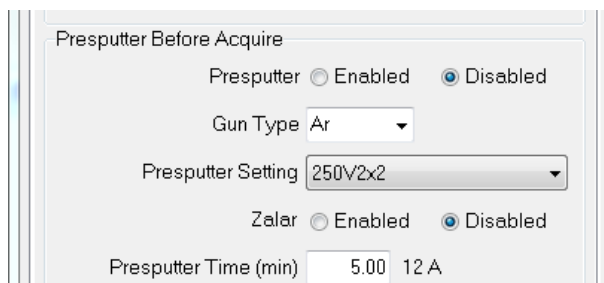

6.9.1. Click 'Enabled' and select the Presputter setting from the drop-down menu for a specific removal rate

Setting	Rate (nm/min)
3kV 2x2	20
2kV 2x2	6.5
1kV 2x2	2.2
500V 2x2	1.3
250V 2x2	0.4

6.9.2. Enter a Presputter time and click 'Close'

6.10 Click on the 'Spe' button the start the survey scan

6.11 Now de-select the box for survey and select all of the other elements for the multiplex analysis

6.12 Click on the 'Spe' button to start the multiplex scan


6.13 All data is immediately available in the PHI Multipak software for processing. See Appendix A for basic data analysis using Multipak, created by Stanford.

7.0 XPS Depth Profile

7.1 Go to the 'XPS' session tab and then the 'Profile' application tab.

7.2 Set up the regions and conditions the same way as in Section 6.

7.3 Click on 'Sputter rate table'

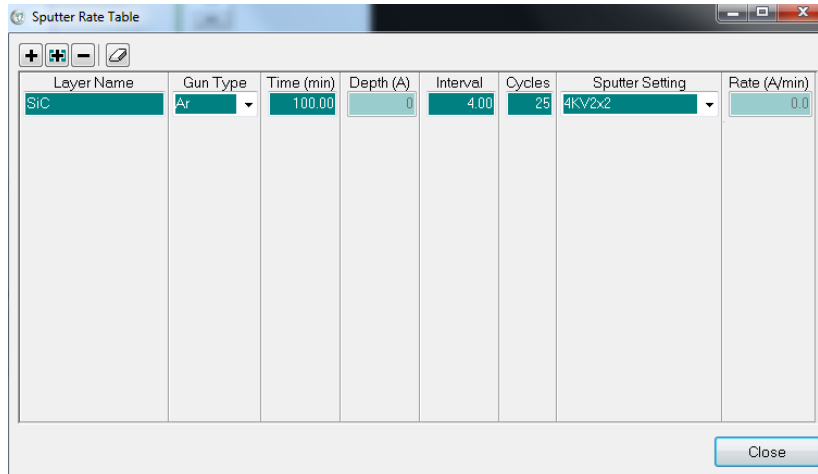

7.4 Select the 'Sputter Setting' from the drop-down menu. The selections and rates are the same as Section 6.

7.5 Enter a time, interval and number of cycles and click on 'Close'.

7.6 Click on the 'Profile' button to begin the scan

8.0 Unloading Sample

8.1 Go to the 'System' session tab and click on the 'Sample transfer' application tab.

8.2 Click on the 'Extract' button.

8.3 After the intro chamber is pumped down, the gate valve will open and a dialog box will appear asking to move the arm in.

8.3.1. Slide the transfer arm all the way until it touches the stop collar and click OK

8.4 The stage will lower the holder on to the transfer arm and another dialog box will appear asking to move the arm out.

8.4.1. Watching through the viewing port, slowly pull the intro rod back completely. V1 should close automatically when the rod is completely withdrawn. Click the OK button in the message dialog to center the stage.

8.4.2. If you see that the sample starts to come off of the arm, then push it all the way back in and try pulling out again.

8.5 The intro chamber will automatically vent. When the lid pops up, use the

metal tweezers and remove the holder.

8.6 While still in the 'System' session tab, click on the 'Vacuum' application tab.

8.7 Select 'Pump intro' from the 'Task' pull-down menu. Then click on 'Process Task'. This will pump down the intro chamber (loadlock chamber).

8.8 Log out of the TUMI

Appendix A: Basic Multipak Analysis (Borrowed from Stanford)

XPS Basic Data Analysis

MultiPak is a MatLab-based program with specific macros written by PHI to handle data reduction of both AES and XPS (ESCA) data.

Once open on the desktop, MultiPak has a "Hide MultiPak", "Show MultiPak" toggle feature that allows you to minimize and restore all the MultiPak windows simultaneously (FIG 1).



FIG 1: Hide/Show shortcut buttons

Spectra are opened in the MultiPak Spectrum window.

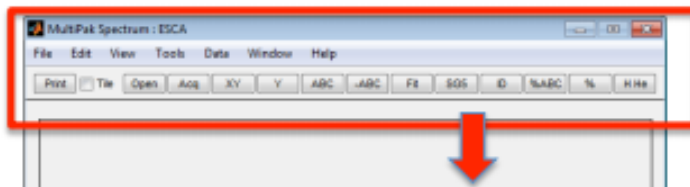


FIG 2: Spectrum Window and shortcut buttons.
* Indicates frequently used.

(1)	(2)	(3)	(4)	(5)	(6)	(7)	(8)	(9)	(10)	(11)	(12)
(1)	Browse and open files										
(2)*	Open last acquisition										
(3)*	Full Scale View										
(4)	View full scale of Y-axis										
(5)	Add annotation										
(6)	Remove annotation										
(7)*	Open Peak Fitting module										
(8)	Applies Savitzky-Golay smoothing to data										
(9)*	Find and Identify Peaks										
(10)*	At% calculation and display										
(11)	Generate at.% data table (text file)										
(12)*	Open Periodic Table										

Task 1: Open and Analyze a Survey Spectrum

1. **Open (1)** or **Acq (2)** to load a spectrum.

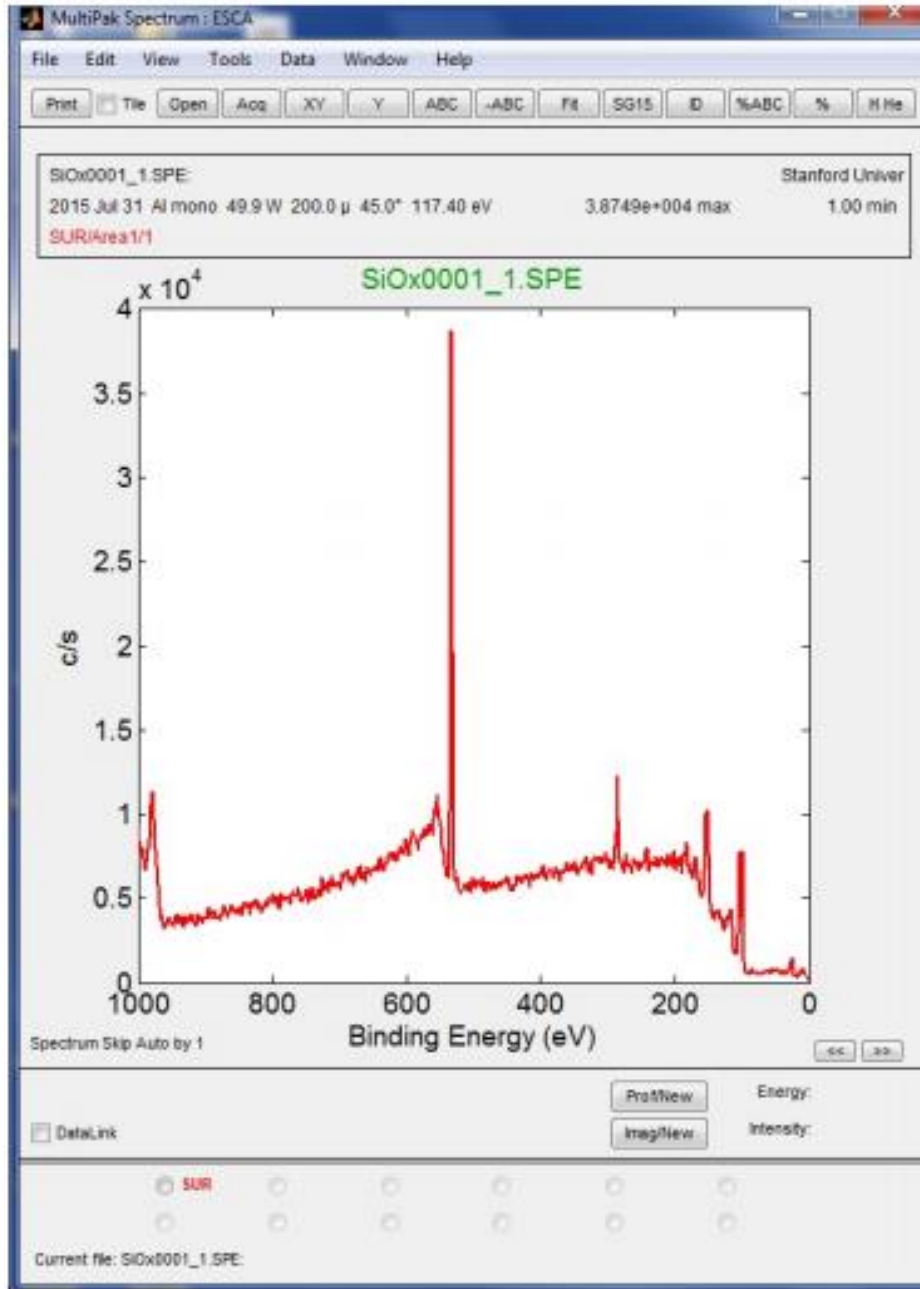


FIG 3. Load a spectrum.

2. Press **ID (9)** to perform Peak Identification. Identified peaks will have labels on the spectrum for display and radio buttons below for atomic composition calculation. See **FIG 4**.

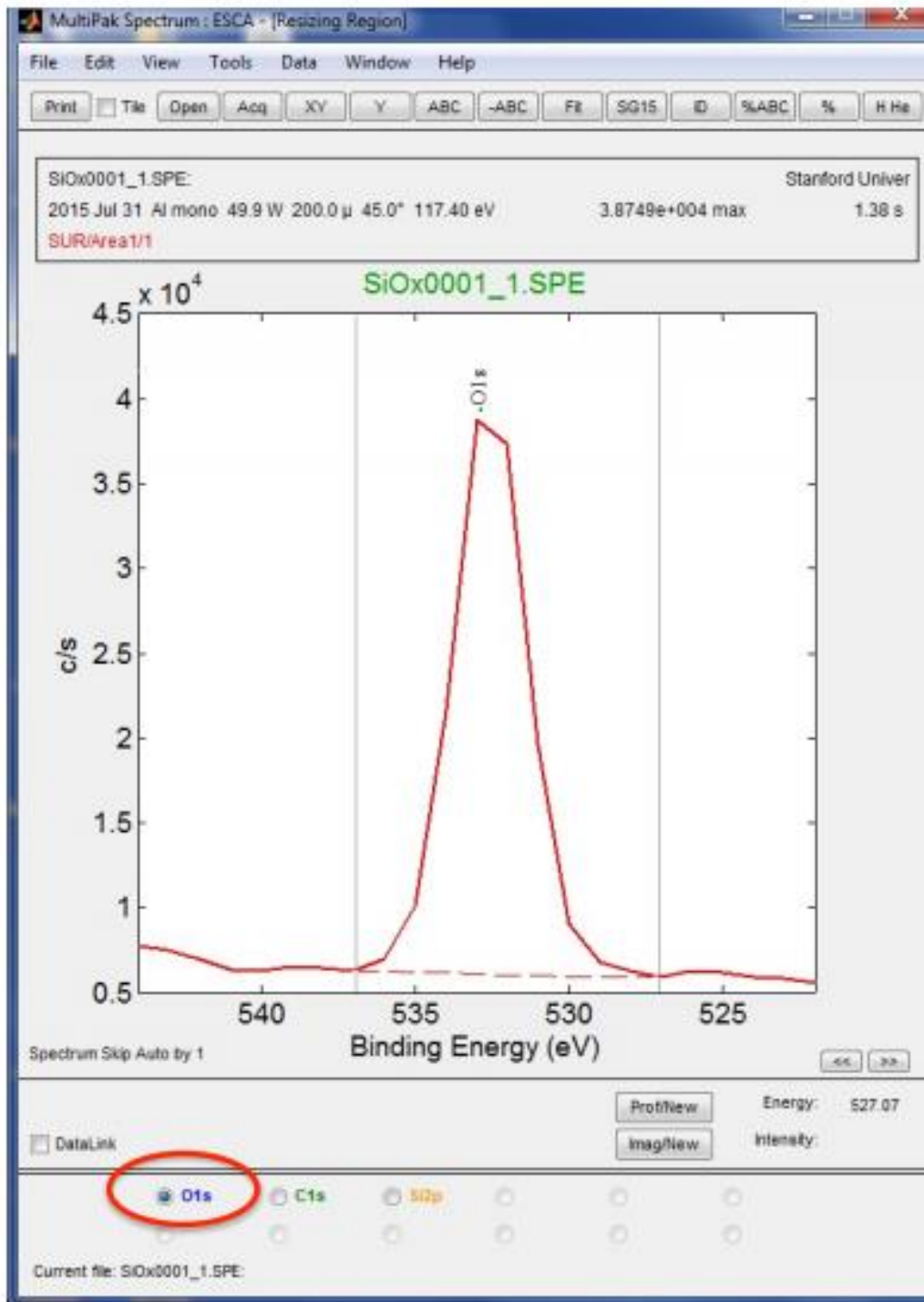


FIG 4. Adjust background.

3. Click on each element transition listed at the bottom of the Spectrum Window to see and adjust the region that will be used to integrate the peak area. Make sure transition label matches the peak(s) (i.e. Ti2p vs. Ti2p3).

4. The identified peaks will also be shown in the Periodic Table window (**FIG 5**). Elements can be added or removed from the spectra and/or atomic percent calculation by clicking on those elements in the Periodic Table.

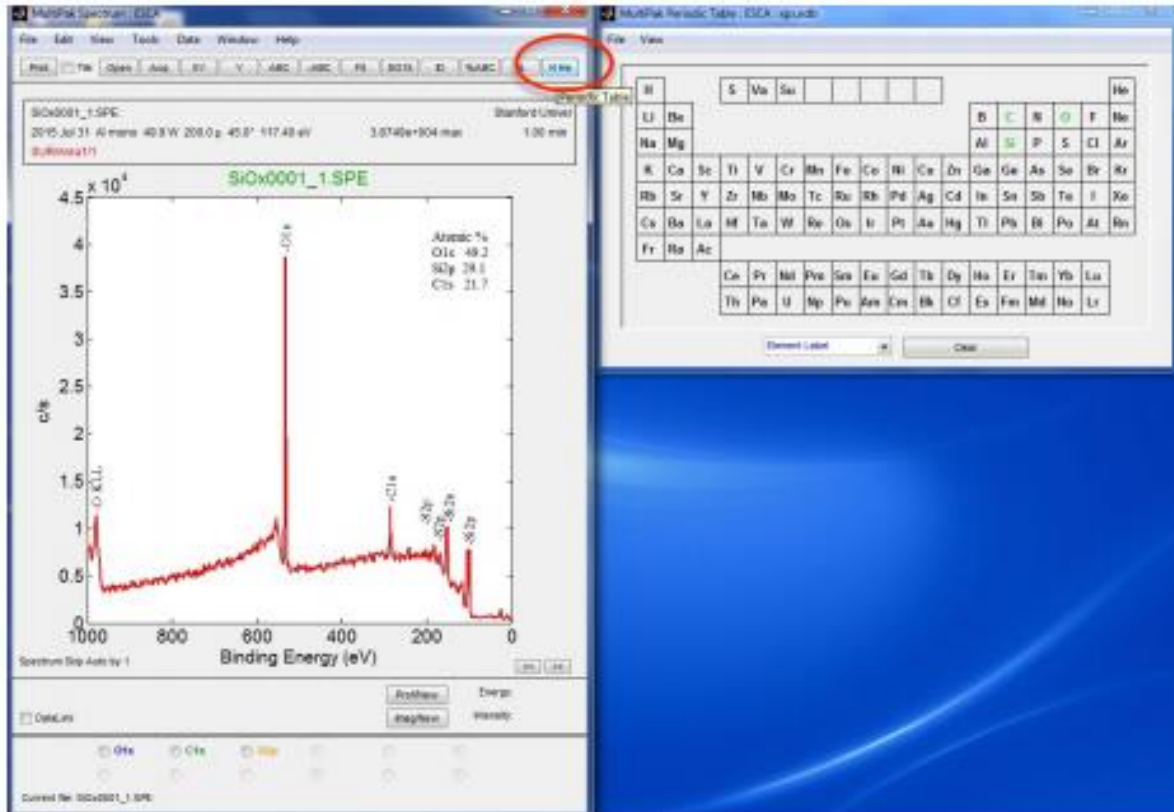


FIG 5: Press H He (12) to open the Periodic Table.

5. For a closer look at a region of the spectrum, zoom by using the left mouse button to click and drag a box. Press **XY (3)** to return to full scale.
6. To actively compare peak positions to known transitions, hold the right mouse button and drag left and right. A green line will mark the binding energy values and transition labels will appear. Release the right mouse button and a desired label will stay on the spectrum.
7. To change the transition of an element used for quantification/labeling:
 - a. Hold SHIFT while selecting the element in the Periodic Table. A Transition dialogue box opens (**FIG 6**).
 - b. In the upper right corner, uncheck the "selected" and "peak label" boxes for the undesired transition.
 - c. Drag down the selector to chose a different transition.
 - d. Recheck the "selected" and "peak label" boxes to make sure a radio button will appear for this transition. When done, press OK.
 - e. Toggle the element off and on in the Periodic Table to apply the change.

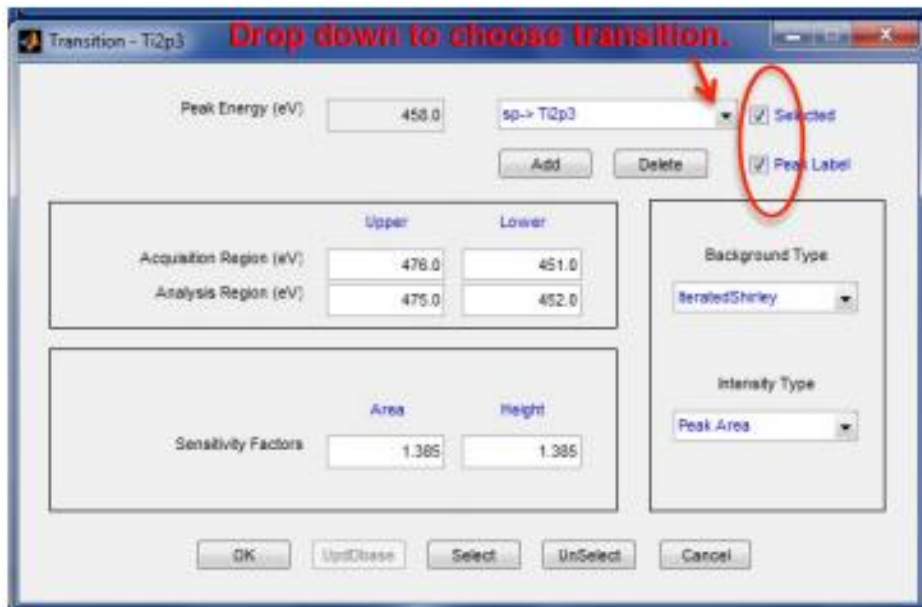


FIG 6. Transition dialogue box.

8. After background fitting is done for all elements listed at the bottom of the Spectrum Window, click the **XY (3)** button to return the graph to Full Scale X and Y.
9. Click the **%ABC (10)** button; the Atomic Concentration table will appear in the upper right corner of the graph.
10. Click the **% (11)** button; the Atomic Concentration table opens in a separate text window that can be saved as a .csv file (**FIG 7**).

```

File Information
-----
File Name: 510x0001_1.SPE
File Type: XPS SPECTRUM
Comment:


Area Information
-----
Area ID: 1
Area Number: 1

Atomic concentration table
-----
Area      Cls      O1s      S12p      RSF      CorrectedRSF
1         0.314    0.733    0.368     49.634   29.10
          36.510   85.530   49.634
          21.71    49.19
  
```

FIG 7. Atomic Concentration Table.

Task 2: Exporting/Saving Data (See FIG 8).

- A. If significant analysis and labeling has been performed go to **File -> Save Current File As** to save the current spectrum as a new .spe file. (Note: Does not save the background subtraction window positions.)
- B. To save current spectrum as a .tiff: **File -> Export To -> tiff.**
- C. To save current spectrum as a .csv: **File -> Export To -> ASCII.**

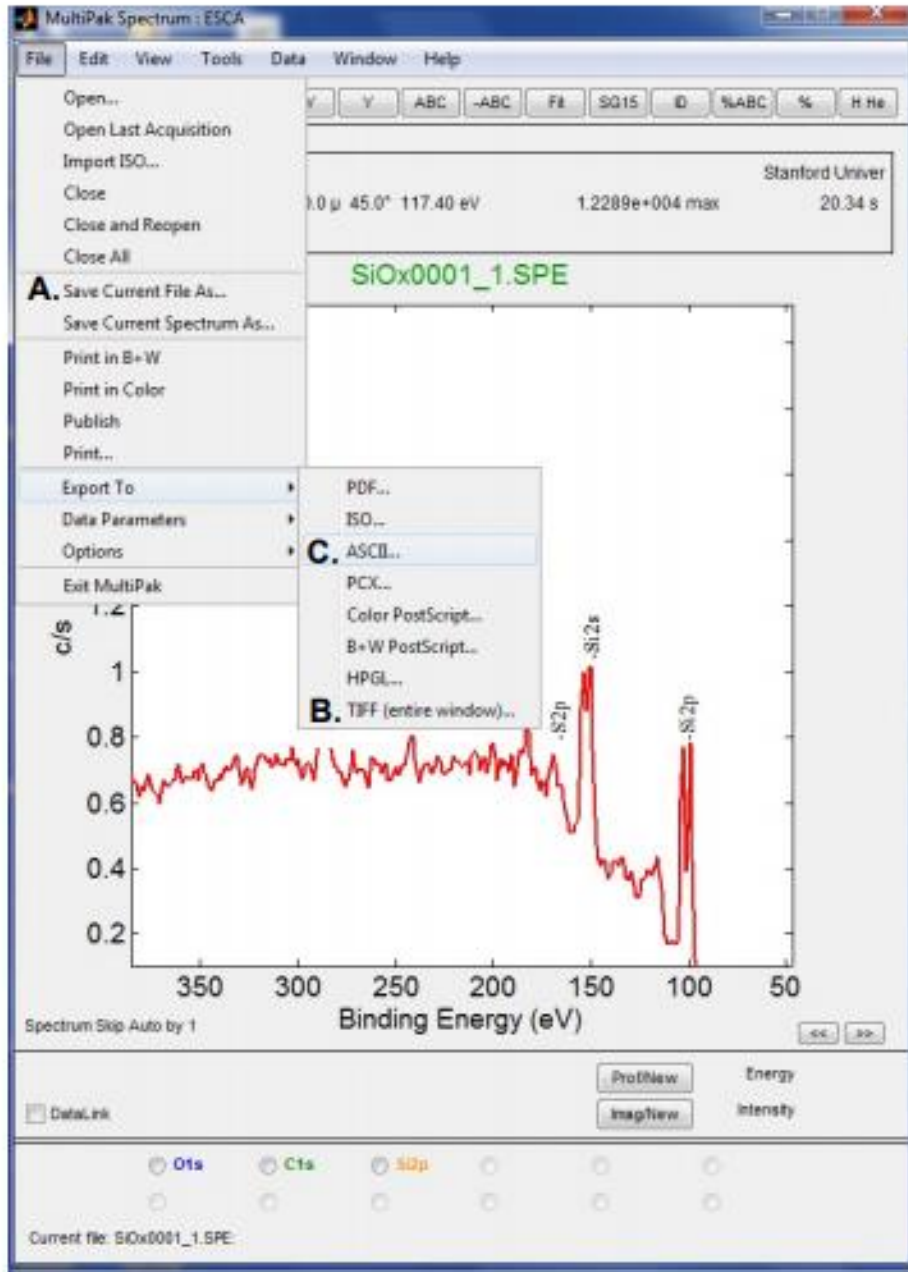**FIG 8. Saving Spectra.**

Task 3: Shifting a Spectrum to Compensate for Charging

Perform this routine on the C1s peak if a spectrum is shifted by a few eV due to charging. If the sample is shifted more than a few eV, use the Neutralizer and Ion Beam Neutralizer in Summit during collection.

1. **Open (1) or Acq (2).**
2. **ID (9)** to Find and Identify Peaks.
3. If the charging has moved the peaks far enough that **ID** does not find the obvious peaks such as C and O, manually click "C" in the Periodic Table Window (**H He (12)**).
4. Click the radio button next for C1s at the bottom of the Spectrum window.
5. In file menu, **Data -> Shift Setup**.
6. A Green line cursor will display in the window. At the bottom of the window it will list PkEnergy and display 284.8eV, the ref. value for C-C/C-H. (See **FIG 9**).

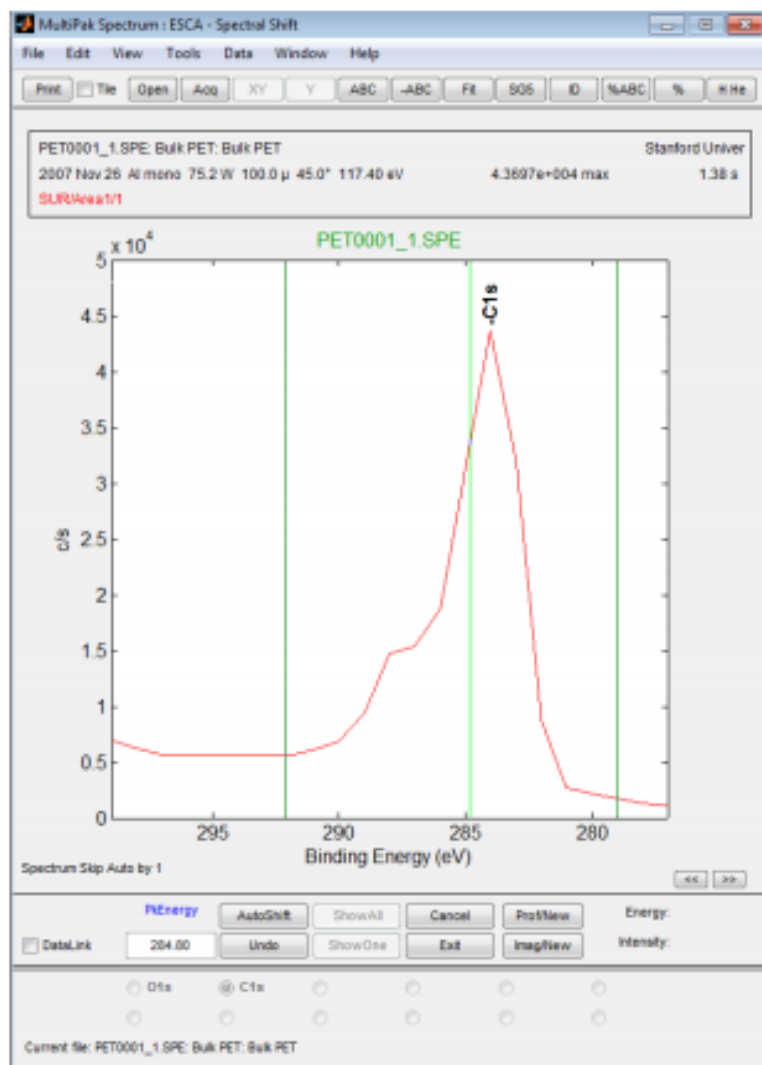


FIG 9. Shift Setup Module in Data.

7. Drag the green cursor to the center of the C-C or C-H peak and the entire spectrum will shift so C1s=284.8eV (FIG 10).

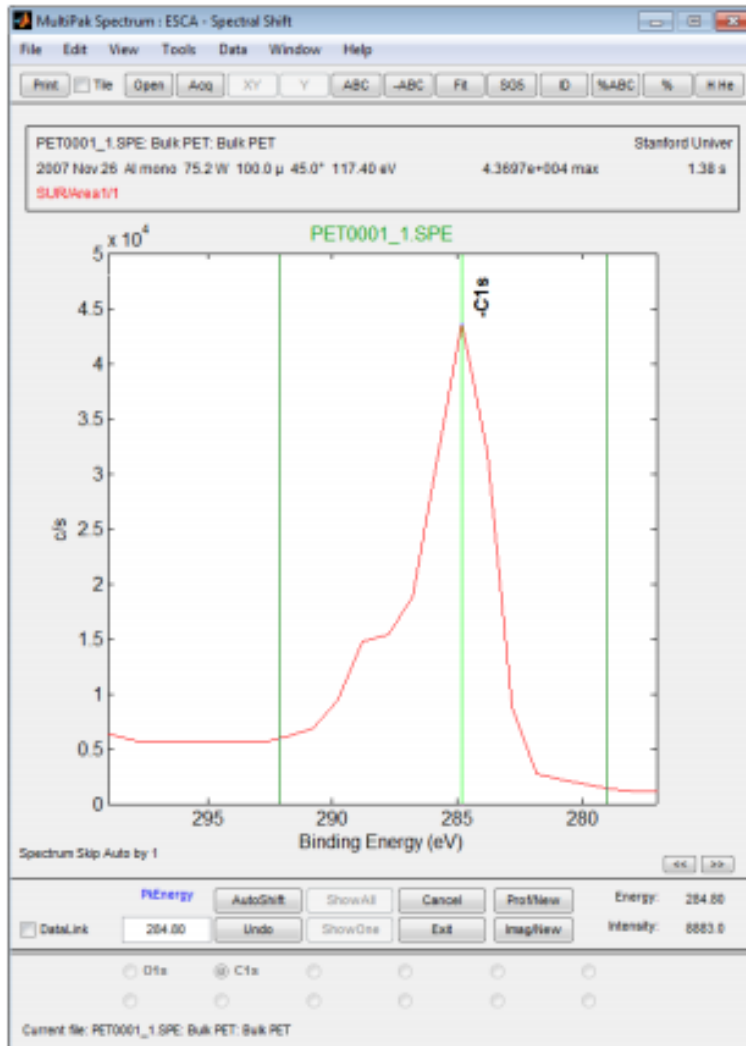


FIG 10. Shifted spectrum. The C1s peak is now centered at the reference value of 284.8eV.

8. If the C1s in some other chemical state such as CF₄ from Teflon, the PkEnergy can be changed to that reference value (for example 292.2eV for Teflon)
9. In the Tools menu, the Chemical State ID feature can be used on a shifted spectrum to search a list of known chemical states.
10. Save the spectrum as a new .spe file since the raw data has been changed (See Task 2).

Task 4: Peak Fitting a High resolution Spectrum

1. **Open (1)** or **Acq (2)** to open a High Resolution spectrum.

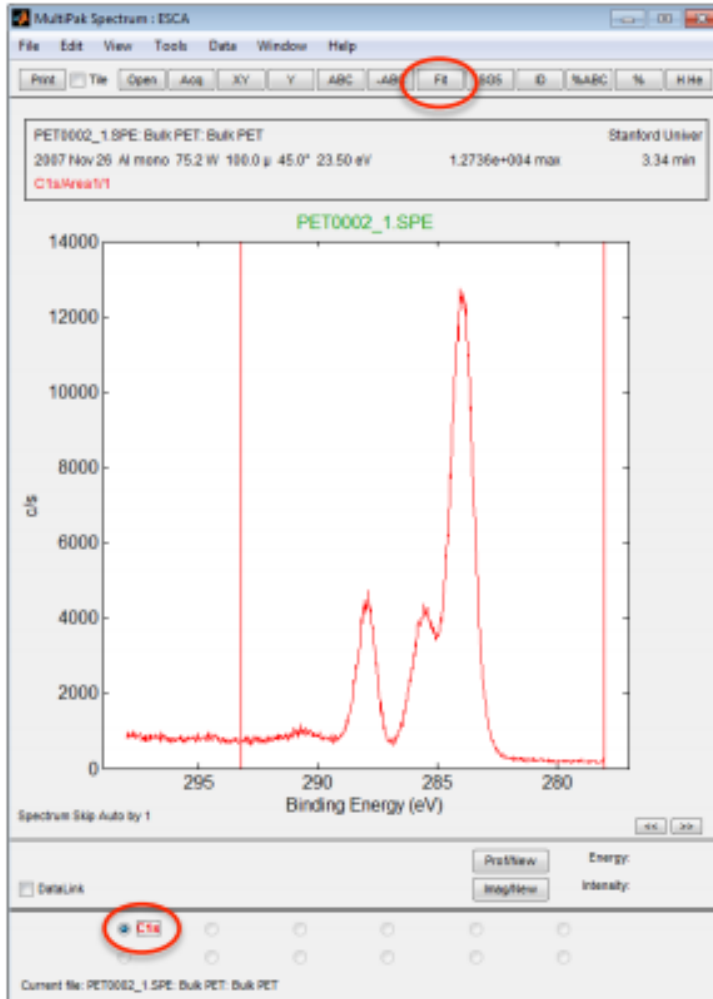


FIG 11. Open a high resolution spectrum. Click FIT.

2. Click the radio button next to the desired element/transition.
3. Click the **FIT (7)** button at the top of the MultiPak Spectrum window. This will open the FIT routine along with an additional Curve Fit Setup window (**FIG 15**). When FIT opens, the previous fit for this element will be applied to the new data. Often the bands (peaks) set up by a previous user will not match your data well.
 - a. See **FIG 12**. Click the **DelAll** button in the bottom of the Spectrum window to clear the fit.
 - b. Adjust background region:
 - i. Uncheck the Setup box in the lower left corner of the Spectrum window.
 - ii. Adjust the left and right had region limits to match the area you wish to fit over.
 - iii. Recheck the Setup box to return to the fit routine.

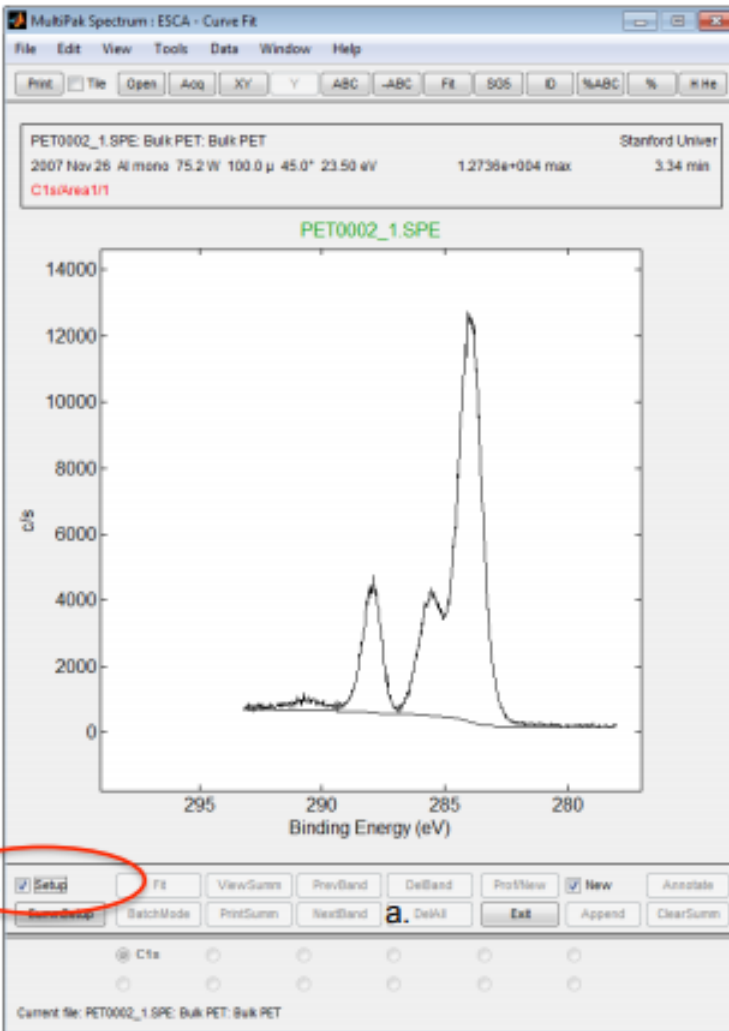


FIG 12. The FIT module after DelAll.

- c. To add a band to the fit, right mouse click in the first peak you want to fit.
 - d. Adjust the band position and width by using the left mouse button on the two square markers on the band (**FIG 13**).
 - e. Repeat (c) and (d) for each peak to be fit.
 - f. Click the **Fit** button in the bottom portion of the Spectrum window.
4. The software will fit the data and display the original data, the bands added for the fit, the fit curve, the background for the fit, and a dotted red line showing the difference between the fit and the original data (Fit Error).

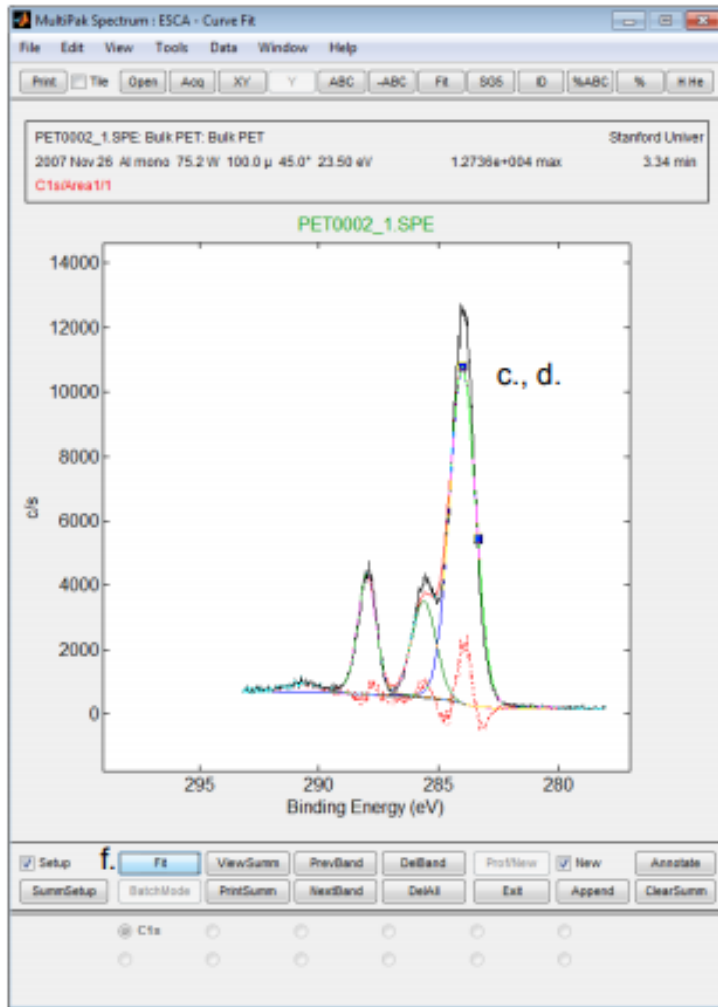


FIG 13. Adding bands to peak fit.

5. The fit results, including χ^2 , are visible in the Curve Fit Setup window (FIG 15).
6. Click the **ViewSum** button at the bottom of the Spectrum window to open a summary (FIG 14) of the fit in a separate text window that can be saved as a .csv file.

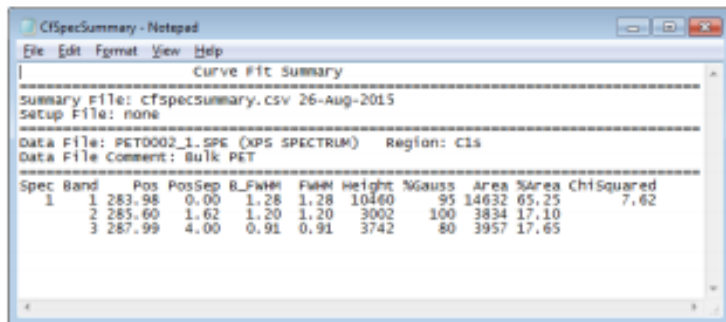


FIG 14. ViewSum button opens summary table.

7. All Fit parameters and some restrictions that can be placed on the fit are shown in the Curve Fit Setup window that opens when you first start the Fit routine.



FIG 15. Curve Fit Setup Window.

8. Saving FIT results:

- a. To save .tiff of curve fit, Spectrum Window -> **File -> Export to-> tiff.**
 - b. To save as individual .spe files to be opened in Spectrum Window:
 - i. In the Curve Fit Setup window (**FIG 15**), click **File->Save Results.**
 - ii. Select your user directory and give the fit result a file name. For example, if the original file name was PET0001_1, then naming the fit result something like PET0001_1Fit is an easy way to keep track of the original data and the fit that goes with it.
 - iii. This will create a series of files that contain each component of the fit. In this case the files are:
 - PET0001_1Fit_1 (First band added)
 - PET0001_1Fit_2 (Second band added)
 - PET0001_1Fit_3 (Third band added)
 - PET0001_1Fit_bg (Background curve for the fit)
 - PET0001_1Fit_err (Error or difference between the fit and original data)
 - PET0001_1Fit_org (Original data curve)
9. Click **Exit** in the bottom of the Spectrum window to exit the fit routine and return to the raw data file.

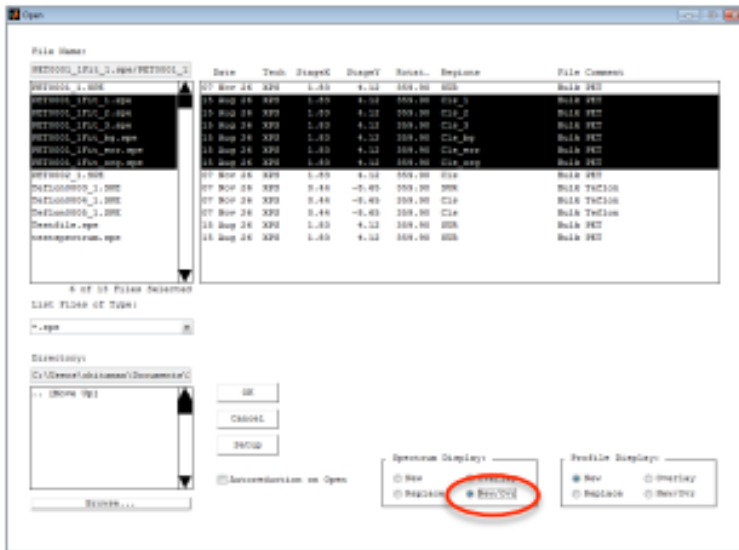


FIG 16. Opening .spe files created by FIT routine.

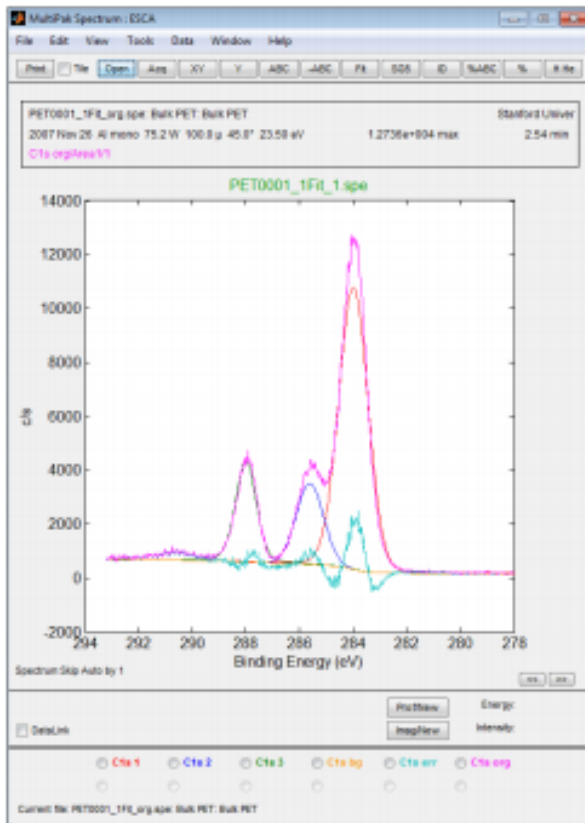


FIG 17. .spe files created in FIT module opened in Spectrum window.

- In Spectrum Window, **Open (1)**. Select all the Files with the PET0001_1Fit_* name and Select **New/Overlay** to overlay all the FIT components in one file. Click **OK (FIG 16)**.
- See Task 2: Exporting/Saving Data to save FIT results as .tiff and .csv files.

Task 5: At.% Composition from FIT Results

Use these steps to consider fitted peaks individually in at.% concentration calculations and/or to separate overlapping peaks from different elements through FIT module.

1. **Open (1)** or **Acq (2)** desired spectrum.
2. Complete Task 3 to shift C1s peak as needed.
3. Complete Task 1 to identify peaks and fit backgrounds without FIT module.
4. Press **Prof/New** at the bottom of the Spectrum window to send intensity data for each element to the Profile window at the right (**FIGs 18, 19**).
5. Perform Task 4 to apply FIT routine to desired peaks.

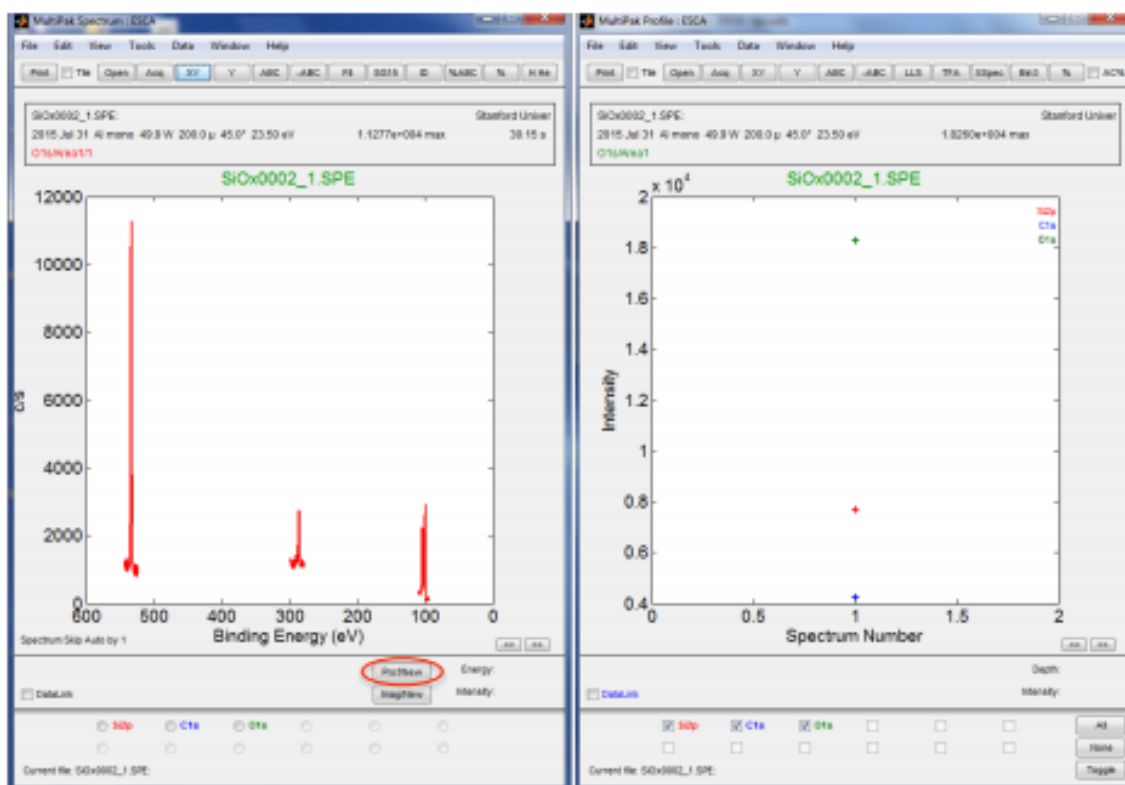


FIG 18: Profile/New Button creates a new Profile window.

6. With the fitted peak in the Spectrum window (FIG 19), uncheck "setup" at the bottom left of the Spectrum window.
7. The **Prof/Update** button will become active (FIG 19). Press **Prof/Update** to send the fitted bands to the Profile window.
8. Note: If a Profile window is open, the **Prof/New** button will change to **Prof/Update**.
9. Repeat Steps 6 & 7 for each FIT result.
10. In the Profile window, at the bottom, uncheck the duplicate elements or bands that should not be included in at.% calculation.
11. Select **AC%** in the upper right corner of the Profile window to display the profile in At.% Composition.

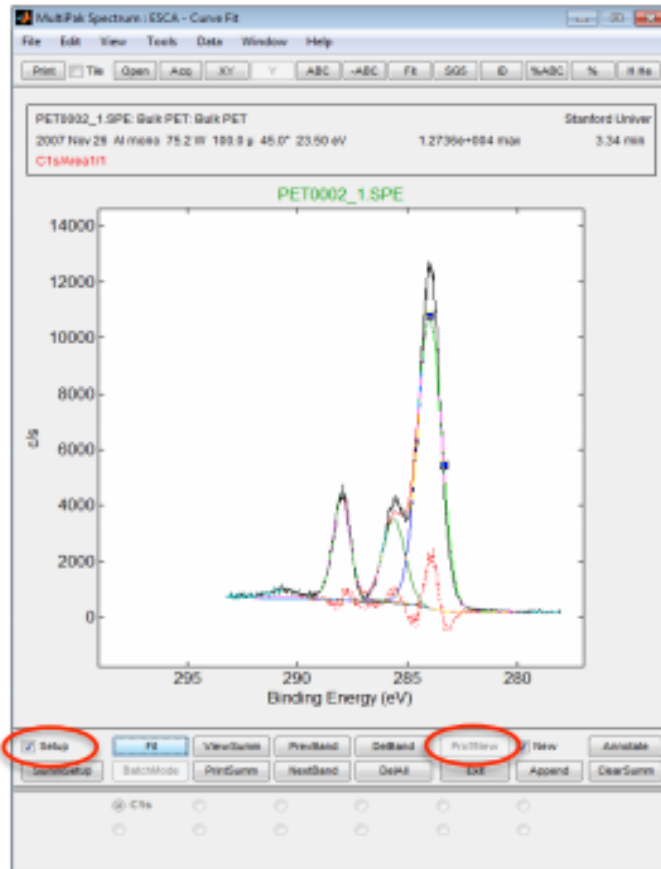


FIG 19. FIT module open in the Spectrum window. Uncheck "setup," press Prof/New or Prof/Update to send bands to Profile window.

12. Press % in the upper right corner of the Profile window to open a text summary of the atomic percent composition results (**FIG 20**).
13. If FIT routine has been used to separate overlapping peaks from different elements, can rename the fitted bands to the correct elements/transitions by holding SHIFT and clicking on the text label in the upper right corner of the profile (**FIG 20**). A Text Properties dialogue box will open. Edit the band to the correct transition, and the atomic percent calculation will use the appropriate sensitivity factor.

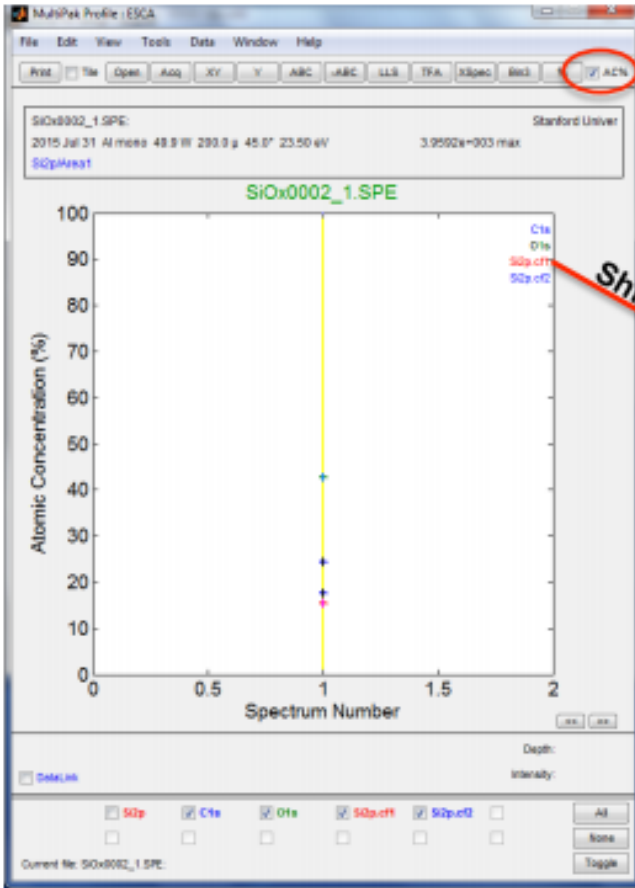


FIG 20. Profile window showing At.% concentration of SiOx sample with bands from Si2p FIT results included (Si2p.cf1, Si2p.cf2). Original Si2p result is unchecked to prevent double counting.

Shift Click

